

Nardis NF-PFP/Biphenyl

Nardis ND-RX

Unison UK-Phenyl









100 x 3 mm

HPLC COLUMN APPLICATION


LC-MS: beta-Carboline alkaloids

β カルボリンアルカロイド類の LC-MS 分析


Nardis NF-PFP/Biphenyl

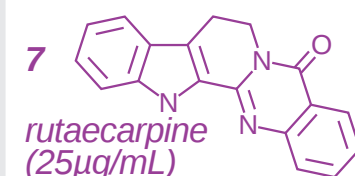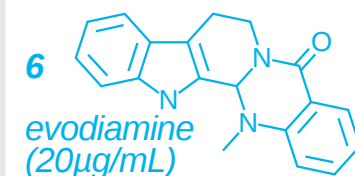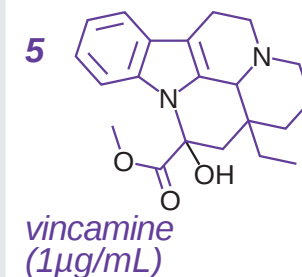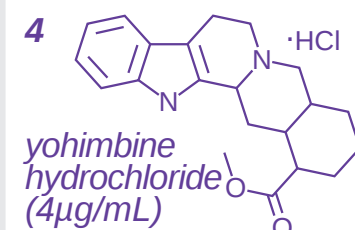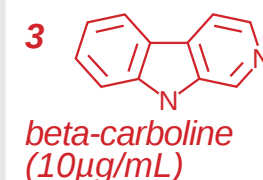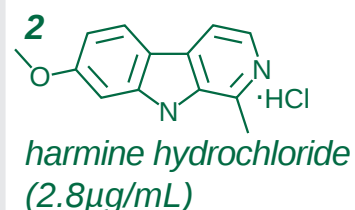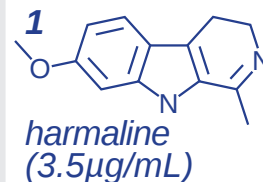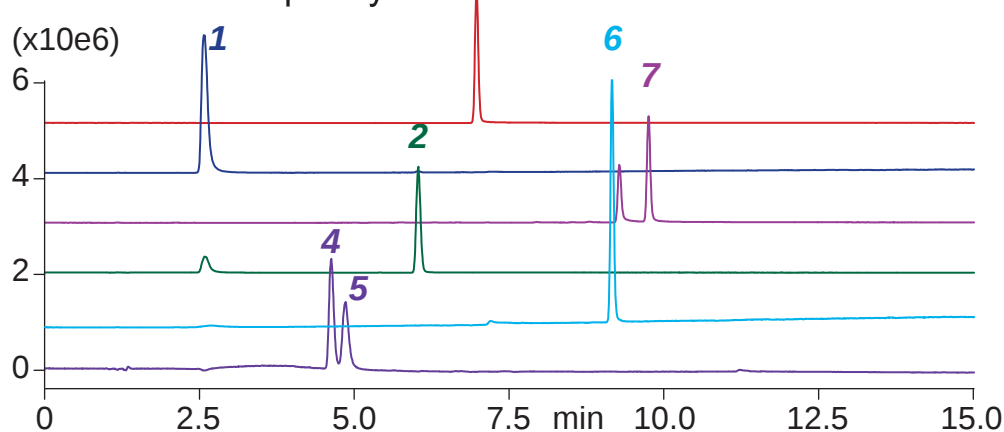

Nardis ND-RX

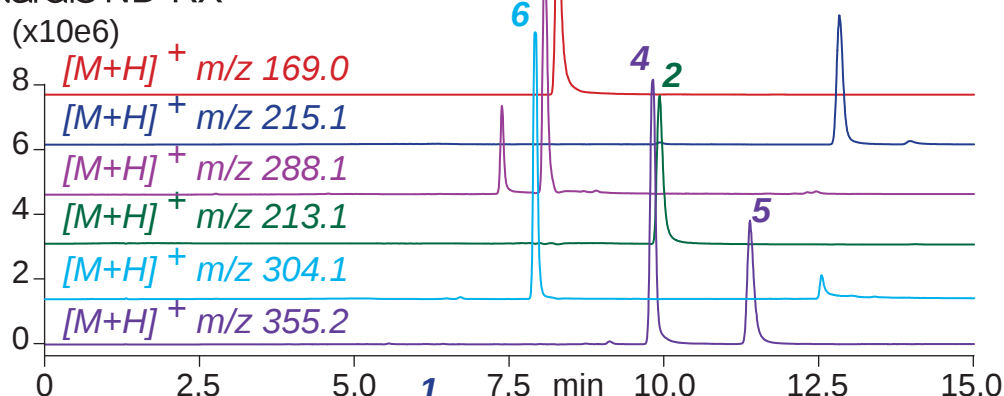

Unison UK-Phenyl

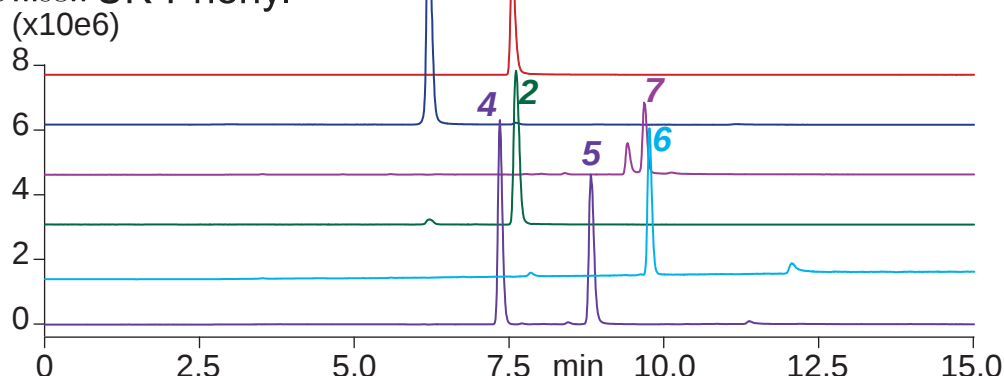

100 x 3 mm

A: acetonitrile / 100mM ammonium acetate = 10 / 90

B: acetonitrile / 100mM ammonium acetate = 80 / 20

0-100%B (0-15min), 100%B (15-17min)

0.4mL/min (7-9MPa), 37°C, Single Quad. MS (ESI)

2μL (1-25μg/mL, ACN / 100mM AcONH₄ = 50 / 50)