

Unison UK-Phenyl


Nardis ND-RX

Nardis NF-PFP/Biphenyl


150 x 4.6 mm

HPLC COLUMN APPLICATION


Isophytol, dihydrophytol(Terpenoid alcohols) イソフィトール、ジヒドロフィトール (テルペノイドアルコール)

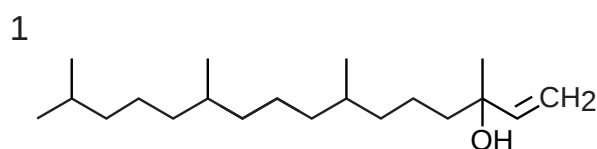

isophytol
(2mg/mL)

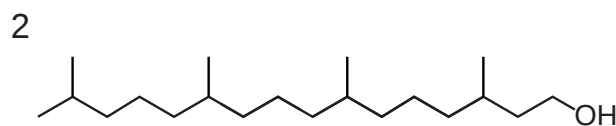

dihydrophytol
(3,7,11,15-tetramethylhexadecan-1-ol,
2mg/mL)

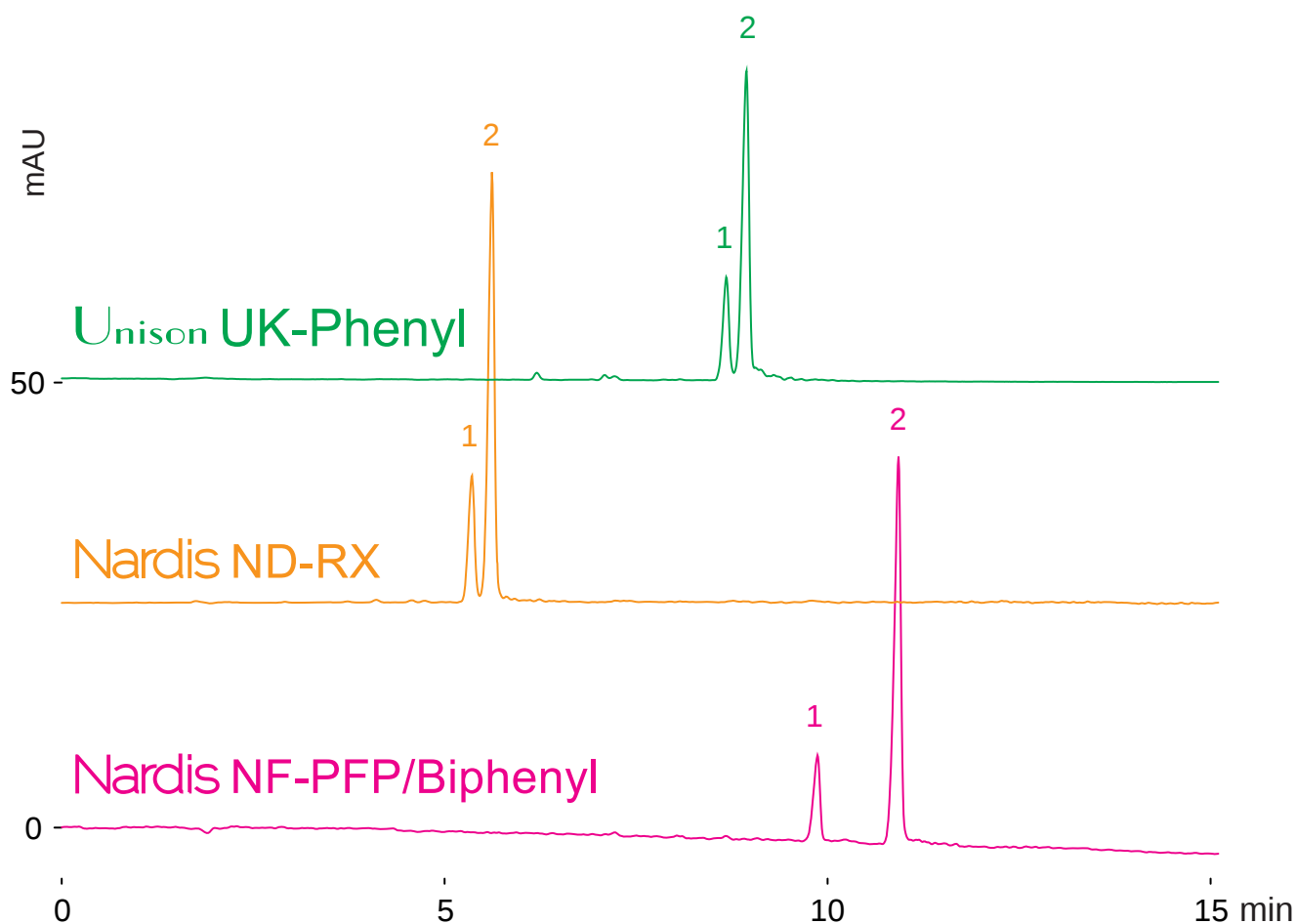

150 x 4.6 mm

A: water / acetonitrile = 50 / 50

B: acetonitrile

0-100%B (0-15min)

1.0mL/min (11-13MPa), 37°C

NQAD (evaporation 35°C, nebulizer 30°C)

10μL (2mg/mL, methanol)