


Scherzo SS-C18

100 x 3 mm

HPLC COLUMN APPLICATION

LC-MS: Lithium ion
リチウムイオンのLC - MS分析

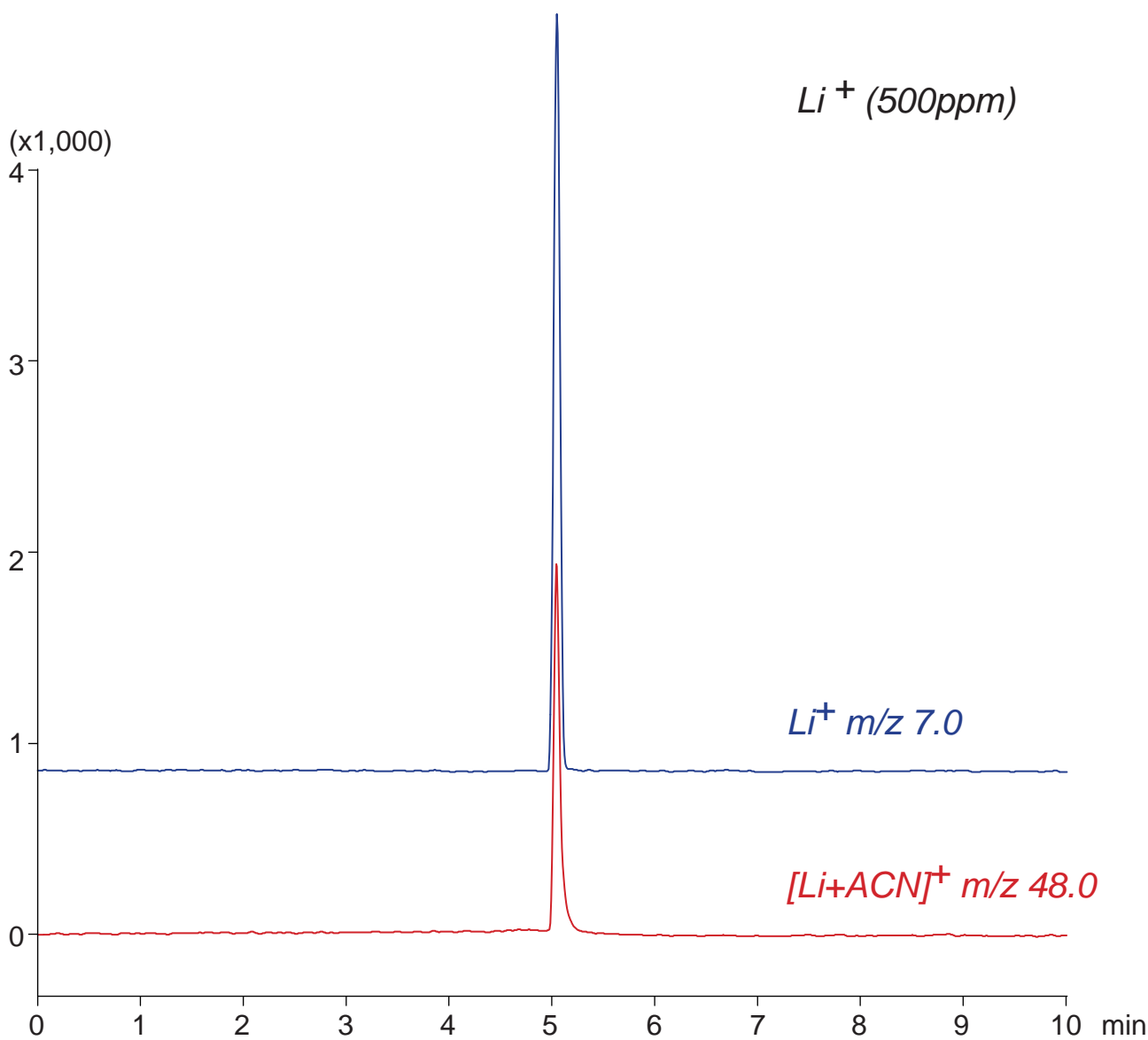

Scherzo SS-C18, 100 x 3 mm

A: water / acetonitrile / formic acid = 10 / 90 / 0.1

B: 100mM ammonium formate / acetonitrile = 80 / 20

0-60%B (0-8min), 100%B (8-10min)

0.4mL/min (8MPa), 37°C, Single Quad. MS (ESI)

5μL (500ppm, water / formic acid = 100 / 1)